



Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

M6399-00	1 x 96 preps
M6399-01	4 x 96 preps
M6399-02	24 x 96 preps

Manual Date: December 2025
Revision Number: v9.0

For Research Use Only

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Table of Contents

Intended Use/Intended User.....	2
Product Description.....	3
Kit Contents/Storage and Stability.....	4
Magnetic Separation Devices and Plasticware.....	4
Preparing Reagents.....	5
Warnings/Safety and Handling Information.....	6
Precautions.....	7
Blood Protocol (250 µL).....	9
Tissue Protocol.....	13
Cultured Cells Protocol.....	18
Saliva Protocol.....	23
Buccal Swabs Protocol.....	27
Mouse Tail Snips Protocol.....	31
Dried Blood Spot Protocol.....	36
Troubleshooting Guide.....	40
Contact Information.....	42
Symbols.....	43
Document Revision History.....	44
Notices and Disclaimers.....	45

Manual Date: December 2025

Revision Number: v9.0



Intended Use

For professional research use.

The Mag-Bind® Blood & Tissue DNA HDQ 96 Kit is intended for isolation and purification of high-quality DNA from whole blood, saliva, mammalian cells and tissue, buccal swabs, mouse tail snips, and dried blood spots.

The Mag-Bind® Blood & Tissue DNA HDQ 96 Kit utilizes magnetic bead-based technology and can be processed either manually or automated on most open-ended liquid handling platforms as well as magnetic processors.

Intended User

The Mag-Bind® Blood & Tissue DNA HDQ 96 Kit is intended for professional use and to be used in a laboratory environment by or under the supervision of professional users, such as laboratory personnel, technicians, researchers, and physicians specifically instructed and trained in molecular biology techniques and familiar with magnetic bead-based purification, either manually or automated.

Product Description

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit offers a versatile method for the isolation of high-quality DNA from a wide variety of samples including fresh or frozen animal cultured cells and tissues, up to 250 µL whole blood, buccal swabs, up to 500 µL saliva, mouse tail snips, and dried blood spots. Mag-Bind® Particles HDQ provide a quick magnetic response time reducing overall processing time. This system combines the reversible nucleic acid-binding properties of Mag-Bind® paramagnetic particles with the time-proven efficiency of Omega Bio-tek's buffer chemistries to provide a fast and convenient method to isolate DNA from a variety of samples. The purification procedure provides high-quality DNA that is suitable for direct use in most downstream applications, such as amplification, next generation sequencing and enzymatic reactions.

If using the Mag-Bind® Blood & Tissue DNA HDQ 96 Kit for the first time, please read this booklet in its entirety to become familiar with the procedures. Samples are lysed in buffer systems that are tailored specifically for each type of starting material. After lysis, samples are mixed with HDQ Binding Buffer and Mag-Bind® Particles HDQ to bind DNA to the magnetic beads. The paramagnetic particles are separated from the lysates by using a magnetic separation device. After a few rapid wash steps to remove trace contaminants, DNA is eluted in Elution Buffer.

Important:

1. If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.
2. Kits include enough reagents for the specified number of preparations plus an additional 10% overage to ensure there is sufficient volume. Please be aware that the actual number of preparations may be lower due to pre-aliquoting of reagents, processing partial plates, and automation platform used etc. Additional reagents are available for purchase separately. Please visit the product page at www.omegabiotek.com or contact your Omega Bio-tek representative for more details and ordering information.

Kit Contents

Product	M6399-00	M6399-01	M6399-02
Preps	1 x 96	4 x 96	24 x 96
Mag-Bind® Particles HDQ	2.2 mL	9 mL	50 mL
AL Buffer	35 mL	125 mL	725 mL
TL Buffer	30 mL	120 mL	600 mL
HDQ Binding Buffer	10 mL	40 mL	200 mL
VHB Buffer	66 mL	230 mL	3 x 440 mL
SPM Buffer	30 mL	150 mL	3 x 150 mL
Elution Buffer	60 mL	250 mL	1000 mL
Proteinase K Solution	2.2 mL	9 mL	50 mL
User Manual	✓	✓	✓

Storage and Stability

All of the Mag-Bind® Blood & Tissue DNA HDQ 96 Kit components are guaranteed for at least 12 months from the date of purchase when stored as follows. Mag-Bind® Particles HDQ should be stored at 2-8°C for long-term use. Proteinase K Solution can be stored at room temperature for up to 12 months. For long-term storage, store Proteinase K Solution at 2-8°C. Store all other components at room temperature and away from bright light. During shipment or storage in cool ambient conditions, precipitates may form in some buffers. Dissolve such deposits by warming the solution at 37°C and gently shaking.

Magnetic Separation Devices and Plasticware

While many brands of magnetic separation devices are compatible with the Mag-Bind® Blood & Tissue DNA HDQ 96 Kit, we recommend using V&P Scientific's Magnetic Plate (Part# VP 771OZM-1) for the binding and wash steps as well as either Alpaqua's Magnum™ EX Magnet Plate (Part# A000380) or Magnum FLX® (Part# A000400) for the elutions steps. These magnetic plates can be used in conjunction with Nunc 2 mL DeepWell™ plates (Part# 278752). This combination provides quick magnetization times, only 1 minute for complete magnetization during wash steps and 5 minutes for lysate clearance steps.

Regardless of the magnetic separation device selected, ensure the device is compatible with the plasticware necessary for this kit.

Preparing Reagents

1. Dilute SPM Buffer with 100% ethanol as follows and store at room temperature.

Kit	100% Ethanol to be Added
M6399-00	70 mL
M6399-01	350 mL
M6399-02	350 mL per bottle

2. Prepare VHB Buffer with 100% ethanol as follows and store at room temperature.

Kit	100% Ethanol to be Added
M6399-00	84 mL
M6399-01	290 mL
M6399-02	560 mL per bottle

3. Prepare HDQ Binding Buffer with 100% isopropanol as follows and store at room temperature.

Kit	100% Isopropanol to be Added
M6399-00	40 mL
M6399-01	160 mL
M6399-02	800 mL

4. Shake or vortex the Mag-Bind® Particles HDQ to fully resuspend the particles before use. The particles must be fully suspended during use to ensure proper binding.

Warnings

This kit is for research use only.

Please read all instructions carefully before using the kit.

Decontaminate and dispose of all potentially infectious materials in accordance with applicable local, state, and national regulations. Please refer to safety data sheets (SDSs) for information on disposal of different components included in this kit.

Safety and Handling Information

All chemicals and biological materials are potentially hazardous.

Biological samples such as plasma, serum, tissues, body fluids, blood etc. are potentially infectious and must be treated as biohazardous materials. Conduct all work in properly equipped facilities following universal precautions and using appropriate personal safety equipment such as disposable gloves, lab coats, safety glasses etc. as required by policies and procedures outlined by your facility. Use appropriate decontaminations and disposal methods in adherence to all applicable local state/provincial, and/or national regulations




Please refer to safety data sheets (SDSs) for information on safe handling, transport, clean-up, and disposal of different components included in this kit. SDSs are made available in PDF format on the product page at www.omegabiotek.com. Discard all waste in accordance with the local safety regulations.

Some of the buffers included in the product contain guanidine-based chaotropic agents, which can form highly reactive compounds when combined with bleach. DO NOT add bleach or acidic solutions to guanidine-containing waste. Please access the SDSs online for detailed information on the reagents.

Where allowed, packaging for non-hazardous buffers, kit boxes, or other packaging materials may be recycled in accordance with local regulations. Reference product labelling or www.omegabiotek.com for guidance.

Precautions

Some of the buffers included in the Mag-Bind® Blood & Tissue DNA HDQ 96 Kit contain guanidine-based chaotropic agents, which can form highly reactive compounds when combined with bleach. **DO NOT add bleach or acidic solutions** to guanidine containing sample-preparation waste. Please access the SDSs online for detailed information on the reagents.

Component	Description
AL Buffer 	Contains: Guanidine hydrochloride. Warning! Causes serious eye irritation. Causes skin irritation. Harmful if swallowed. Do not eat, drink or smoke when using this product. Wash all exposed external body areas thoroughly after handling. Wear protective gloves, protective clothing, eye protection, and face protection. IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Get medical advice/attention if eye irritation persists. Take off contaminated clothing and wash before reuse. ON SKIN: Wash with plenty of water and soap. Get medical advice/attention if skin irritation or rash occurs. SWALLOWED: Rinse mouth. Call a poison center or doctor/physician if you feel unwell.
TL Buffer 	Contains: Anionic detergent. Warning! Causes serious eye irritation. May cause an allergic skin reaction. Avoid breathing mist/vapors/spray. Contaminated work clothing should not be allowed out of the workplace. Wear protective gloves/protective clothing/eye protection/face protection. IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Get medical advice/attention if eye irritation persists. ON SKIN: Wash with plenty of water and soap. Get medical advice/attention if skin irritation or rash occurs. Wash contaminated clothing before reuse.
Proteinase K Solution 	Contains: Proteinase K. Danger! Causes mild skin irritation. May cause allergy or asthma symptoms or breathing difficulties if inhaled. Avoid breathing dust/fume/gas/mist/vapors/spray. Wear protective gloves/protective clothing/eye protection/face protection. Wear respiratory protection. If exposed or concerned: Call a poison center or doctor/physician. Remove victim to fresh air and keep at rest in a position comfortable for breathing.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Blood Protocol (250 µL)

The procedure below has been optimized for use with 250 µL FRESH or FROZEN blood samples. Buffy coat can also be used.

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Reagents to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Heat block, incubator, or water bath capable of 70°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plate (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- Optional: RNase A (10 mg/mL)
- Optional: PBS

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
 - Set heat block, incubator, or water bath to 70°C.
1. Prepare a mastermix of AL Buffer and Proteinase K Solution only for samples to be extracted according to the table below:

Component	Amount per Prep	Total Amount per 96-well Plate
AL Buffer	290 µL	30.6 mL*
Proteinase K Solution	20 µL	2.1 mL*

* 10% excess volume has been calculated for a 96-well plate.

Important:

- Only prepare as much AL Buffer/Proteinase K Solution master mix that will be used within 4 hours of preparation.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

- Mix the AL Buffer/Proteinase K Solution master mix by inverting carefully.
 - Remove any bubbles from the AL Buffer/Proteinase K Solution master mix if using with automated systems.
2. Add 250 μ L blood sample to a 2 mL 96-well deep-well plate (not provided). If the volume of blood is less than 250 μ L, bring the volume up to 250 μ L with PBS (not provided) or Elution Buffer.

Note: For buffy coat samples, we recommend starting with 100 μ L and bringing the volume up to 250 μ L with PBS (not provided) or Elution Buffer.

3. Add 310 μ L AL Buffer/Proteinase K Solution master mix to each sample. Vortex or pipet up and down 20 times to mix. Proper mixing is crucial for good yield.

Note: For automated protocols tip mixing yields best results and is recommended.

4. Incubate for 10 minutes at 70°C.

Optional: Add 5 μ L RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

5. Add 400 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes

6. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
7. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
8. Remove the plate from the magnetic separation device.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

9. Add 600 μ L VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

10. Vortex for 2 minutes.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

11. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

12. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

13. Remove the plate from the magnetic separation device.

14. Repeat Steps 9-13 for a second VHB Buffer step.

15. Add 600 μ L SPM Buffer.

Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

16. Vortex for 2 minutes.

17. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

18. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
19. Select one of the following ethanol removal steps:
 - A. Leave the plate on the magnetic separation device. Add 500 µL nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 20.

OR

- B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 20.
20. Remove the plate from the magnetic separation device.
21. Add 50-200 µL Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.

22. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.

23. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
24. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).
25. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Tissue Protocol

This method allows genomic DNA isolation from up to 10 mg tissue. Yields will vary depending on the source.

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Equipment to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Centrifuge with swing-bucket rotor capable of 4,000g
- Centrifuge adapter for 96-well plates
- Shaking water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plates (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- Recommended: 1M Dithiothreitol (DTT)
- Optional: RNase A (10 mg/mL)
- Optional: Heat block, incubator, or water bath capable of 70°C
- Optional: Liquid nitrogen and mortar and pestle

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
- Set water bath to 55°C.
- Optional: Set water bath, incubator, or heat block 70°C.
- Recommended: Add 40 µL 1M DTT per 1 mL TL Buffer before use.

OPTIONAL: Although mechanical homogenization of tissue is not necessary, pulverizing the samples in liquid nitrogen will improve lysis and reduce incubation time. Once the liquid nitrogen has evaporated, transfer the powdered tissue to a clean 96-well deep-well plate (not provided). Add 250 µL TL Buffer and proceed to Step 3 on the next page.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

1. Mince up to 10 mg tissue and transfer to a 96-well deep-well plate (not provided).

Note: Cutting the tissue into small pieces can speed up lysis.

2. Add 250 μ L TL Buffer.

Optional: For lysis of hair or other tough-to-lyse tissues, a mastermix of TL Buffer and DTT is recommended.

- Dilute DTT to a final concentration of 40 mM in TL Buffer.
- Add 40 μ L 1M DTT per 1 mL TL Buffer before use.
- Only prepare as much TL Buffer/DTT mastermix that will be used immediately.

3. Add 20 μ L Proteinase K Solution. Vortex to mix.

4. Incubate at 55°C in a shaking water bath.

Note: If a shaking water bath is not available, vortex the sample every 20-30 minutes. Lysis time depends on amount and type of tissue, but is usually under 3 hours. The lysis can proceed overnight.

Optional: Add 5 μ L RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

5. Centrifuge at maximum speed ($\geq 4,000g$) for 5 minutes to pellet undigested tissue debris.
6. Carefully transfer 200 μ L of the supernatant to a new 96-well deep-well plate without disturbing the undigested pellet.
7. Add 230 μ L AL Buffer to each sample. Vortex for 10 minutes to mix. Proper mixing is crucial for good yield.

Note:

- For automated protocols tip mixing yields best results and is recommended.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

8. Add 320 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ to each sample. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

9. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
10. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
11. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
12. Add 600 μ L VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

13. Vortex for 2 minutes to mix.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

14. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
15. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

16. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
17. Repeat Steps 12-16 for a second VHB Buffer step.
18. Add 600 µL SPM Buffer.

Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
19. Vortex for 2 minutes to mix.
20. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
21. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
22. Select one of the following ethanol removal steps:
 - A. Leave the plate on the magnetic separation device. Add 500 µL nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 23.

OR

 - B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 23.
23. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

24. Add 100-200 μ L Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.

25. Vortex for 5 minutes to mix.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.

26. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

27. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).

28. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Cultured Cells Protocol

This protocol is designed for rapid isolation of up to 25 µg genomic DNA from up to 5 x 10⁶ cultured cells.

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Equipment to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Centrifuge with swing-bucket rotor capable of 4,000g
- Shaking water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plates (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- Cold PBS (4°C)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- Optional: RNase A (10 mg/mL)
- Optional: Heat block, incubator, or water bath capable of 70°C
- Optional: Trypsin and cell scraper

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
- Set shaking water bath to 55°C.
- Optional: Set water bath, incubator, or heat block 70°C.

1. Prepare the cell suspension.

- 1a. Frozen cell samples should be thawed before starting this protocol. Pellet cells by centrifugation. Wash the cells with cold PBS (4°C) and resuspend cells in 180 µL cold PBS (not provided). Proceed with Step 2 of this protocol.
- 1b. For cells grown in suspension, pellet 5 x 10⁶ cells at 1,200g in a centrifuge

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

tube. Discard the supernatant, wash the cells once with cold PBS (4°C), and resuspend cells in 180 µL cold PBS (not provided). Proceed with Step 2 of this protocol.

- 1c. For cells grown in a monolayer, harvest the cells by either using a trypsin treatment or cell scraper. Wash cells twice in cold PBS (4°C) and resuspend the cells with 180 µL cold PBS. Proceed with Step 2 of this protocol.
2. Prepare a mastermix of AL Buffer and Proteinase K Solution only for samples to be extracted according to the table below:

Component	Amount per Prep	Total Amount per 96-well Plate
AL Buffer	230 µL	24.3 mL*
Proteinase K Solution	20 µL	2.1 mL*

* 10% excess volume has been calculated for 96-well plate.

Important:

- Only prepare as much AL Buffer/Proteinase K Solution master mix that will be used within 4 hours of preparation.
 - Mix the AL Buffer/Proteinase K Solution master mix by inverting carefully.
 - Remove any bubbles from the AL Buffer/Proteinase K Solution master mix if using with automated systems.
3. Add 250 µL AL Buffer/Proteinase K Solution mastermix to each sample. Vortex for 10 minutes to mix. Proper mixing is crucial for good yield.

Note:

- For automated protocols tip mixing yields best results and is recommended.
 - If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.
4. Incubate for 10 minutes at 55°C in a shaking water bath.

Note: If a shaking water bath is not available, vortex the samples every 2-3 minutes.

5. Transfer the samples to a 96-well deep-well plate (not provided).

Optional: Add 5 µL RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

6. Add 320 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ. Vortex for 10 minutes.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
 - If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.
7. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
 8. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
 9. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
 10. Add 600 μ L VHB Buffer.
Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
 11. Vortex for 2 minutes.
Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.
 12. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
 13. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
 14. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

15. Repeat Steps 10-14 for a second VHB Buffer step.
16. Add 600 μ L SPM Buffer.
Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
17. Vortex for 2 minutes to mix.
18. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
19. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
20. Select one of the following ethanol removal steps:

- A. Leave the plate on the magnetic separation device. Add 500 μ L nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 21.

OR

- B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 21.
21. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
22. Add 50-200 μ L Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

23. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.

24. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

25. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).

26. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Saliva Protocol

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Equipment to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Shaking water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plates (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- Optional: RNase A (10mg/mL)
- Optional: Heat block, incubator, or water bath capable of 70°C

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
- Set shaking water bath to 55°C.
- Optional: Set water bath, incubator, or heat block 70°C.

1. Centrifuge the saliva tube at 2,000g for 5 minutes.
2. Transfer 500 µL stabilized saliva samples (e.g. DNA Genotek Oragene®, Mawi iSWAB™, Biomatrix® DNAGard® Saliva) to a 96-well deep-well plate (not provided).
3. Prepare a mastermix of AL Buffer and Proteinase K Solution only for samples to be extracted according to the table below:

Component	Amount per Prep	Total Amount per 96-well Plate
AL Buffer	200 µL	21.12 mL*
Proteinase K Solution	20 µL	2.1 mL*

* 10% excess volume has been calculated for 96-well plate.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Important:

- Only prepare as much AL Buffer/Proteinase K Solution master mix that will be used within 4 hours of preparation.
- Mix the AL Buffer/Proteinase K Solution master mix by inverting carefully.
- Remove any bubbles from the AL Buffer/Proteinase K Solution master mix if using with automated systems.

4. Add 220 μ L AL Buffer/Proteinase K Solution to each sample. Vortex for 10 minutes to mix. Proper mixing is crucial for good yield.

Note:

- For automated protocols tip mixing yields best results and is recommended.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

5. Incubate for 10 minutes at 55°C in a shaking water bath.

Note: If a shaking water bath is not available, vortex the plate every 2-3 minutes. If DNA Genotek Oragene® tube was used and incubation step was already performed, skip to Step 6.

Optional: Add 5 μ L RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

6. Add 400 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

7. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

8. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

9. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

10. Add 600 μ L VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

11. Vortex for 2 minutes.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

12. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

13. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

14. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.

15. Repeat Steps 10-14 for a second VHB Buffer step.

16. Add 600 μ L SPM Buffer.

Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

17. Vortex for 2 minutes.

18. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

19. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
20. Select one of the following ethanol removal steps:
 - A. Leave the plate on the magnetic separation device. Add 500 µL nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 21.

OR

- B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 21.
21. Add 100-200 µL Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.
22. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.
23. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
24. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).
25. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Buccal Swabs Protocol

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Equipment to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Centrifuge with swing-bucket rotor capable of 4,000g
- Centrifuge adapter for 96-well plates
- Shaking water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plates (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Optional: RNase A (10 mg/mL)
- Optional: Nuclease-free water
- Optional: Heat block, incubator, or water bath capable of 70°C

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
 - Set shaking water bath to 55°C.
 - Optional: Set water bath, incubator, or heat block 70°C.
1. Cut off the buccal brush or swab head and place each swab into a well of a 96-well deep-well plate (not provided).

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

2. Prepare a mastermix of AL Buffer, Proteinase K Solution, and Elution Buffer only for samples to be extracted according to the table below:

Component	Amount per Prep	Total Amount per 96-well Plate
AL Buffer	290 μ L	30.6 mL*
Proteinase K Solution	20 μ L	2.1 mL*
Elution Buffer	250 μ L	26.4 mL

* 10% excess volume has been calculated for a 96-well plate.

Important:

- Only prepare as much AL Buffer/Proteinase K Solution master mix that will be used within 4 hours of preparation.
 - Mix the AL Buffer/Proteinase K Solution master mix by inverting carefully.
 - Remove any bubbles from the AL Buffer/Proteinase K Solution master mix if using with automated systems.
3. Add 560 μ L AL Buffer/Proteinase K Solution/Elution Buffer mastermix to each sample. Vortex or pipet up and down 20 times to mix.
- Note:** For automated protocols tip mixing yields best results and is recommended.
4. Incubate for 10 minutes at 55°C in a shaking water bath.
- Note:** If a shaking water bath is not available, vortex the plate every 2-3 minutes.
5. Centrifuge at 3,000g for 2 minutes.
6. Transfer 500 μ L lysate into a new 96-well deep-well plate. Do not transfer the swabs to the new plate.

Optional: Add 5 μ L RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

7. Add 350 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

8. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
9. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
10. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
11. Add 600 µL VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

12. Vortex for 2 minutes.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

13. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
14. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
15. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
16. Repeat Steps 11-15 for a second VHB Buffer step.
17. Add 600 µL SPM Buffer.

Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

18. Vortex for 2 minutes.

19. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

20. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

21. Leave the plate on the magnetic separation device for 10 minutes to air dry the Mag-Bind® Particles HDQ. Remove any residual liquid from the wells.

Note: All liquid must be aspirated at this step. It is helpful to remove all liquid from the well then wait one minute and remove any residual liquid from the well.

22. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.

23. Add 100-200 µL Elution Buffer or nuclease-free water (not provided) to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.

24. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.

25. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

26. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).

27. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Mouse Tail Snips Protocol

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Equipment to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Centrifuge with swing-bucket rotor capable of 4,000g
- Centrifuge adapter for 96-well plates
- Shaking water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plates (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- Recommended: 1M Dithiothreitol (DTT)
- Optional: RNase A (10 mg/mL)
- Optional: Heat block, incubator, or water bath capable of 70°C

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
 - Set water bath to 55°C.
 - Recommended: Add 40 µL 1M DTT per 1 mL TL Buffer.
1. Snip a 2-5 mm piece of mouse tail, cut into several pieces, and transfer the pieces to a 96-well deep-well plate (not provided).

Note: Follow all regulations regarding the safe and humane treatment of animals. Mice should not be older than 6 weeks since lysis will be more difficult in older animals resulting in suboptimal DNA yields. If possible, obtain tail biopsies at 2-4 weeks and freeze samples at -70°C until DNA is extracted.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

2. Add 250 μ L TL Buffer.

Optional: For lysis of hair or other tough-to-lyse tissues, a mastermix of TL Buffer and DTT is recommended.

- Dilute DTT to a final concentration of 40 mM in TL Buffer.
- Add 40 μ L 1M DTT per 1 mL TL Buffer before use.
- Only prepare as much TL Buffer/DTT mastermix that will be used immediately.

3. Add 20 μ L Proteinase K Solution to each sample. Vortex to mix.

4. Incubate at 55°C in a shaking water bath for 1-4 hours or until lysis is complete.

Note:

- If a shaking water bath is not available, vortex the samples vigorously every 20-30 minutes. Incomplete lysis may significantly reduce DNA yields. Incubation time for complete tail lysis is dependent on length of tail snip and age of animal, e.g. a 5 mm tail piece from a 2 week old mouse typically will lyse in 2 hours.
- For older animals, an overnight incubation may improve yields. Note that bone and hair will not lyse. Adding DTT to TL Buffer at a final concentration of 40 mM can help with tissue lysis.

5. Centrifuge at maximum speed ($\geq 4,000g$) for 5 minutes to pellet undigested tissue debris.

6. Carefully transfer 200 μ L supernatant to a new 96-well deep-well plate without disturbing the undigested pellet.

Optional: Add 5 μ L RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

7. Add 230 μ L AL Buffer. Vortex for 10 minutes to mix. Proper mixing is crucial for good yield.

Note:

- For automated protocols tip mixing yields best results and is recommended.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

8. Add 320 μ L HDQ Binding Buffer and 20 μ L Mag-Bind® Particles HDQ. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes.

9. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
10. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
11. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
12. Add 600 μ L VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

13. Vortex for 2 minutes.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

14. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
15. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

16. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
17. Repeat Steps 12-16 for a second VHB Buffer step.
18. Add 600 μ L SPM Buffer.

Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
19. Vortex for 2 minutes.
20. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
21. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
22. Select one of the following ethanol removal steps:
 - A. Leave the plate on the magnetic separation device. Add 500 μ L nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 23.

OR

 - B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 23.
23. Remove the plate containing the Mag-Bind® Particles HDQ from the magnetic separation device.
24. Add 100-200 μ L Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

25. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.

26. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.

27. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).

28. Store DNA at -20°C.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

Dried Blood Spot Protocol

Important: If automating this procedure on a liquid handler or a magnetic processor, please contact your Omega Bio-tek representative for instrument-specific instructions. It is the responsibility of the user to validate any automated method for any particular use.

Materials and Reagents to be Supplied by User:

- Magnetic separation device - Recommend dual magnet approach:
 - Bind/wash steps: V&P Scientific, Part#VP 771OZM-1
 - Elution step: Either Alpaqua Magnum™ EX, Part#A000380 or Magnum FLX® Part#A000400
- Vortexer
- Heat block, incubator, or water bath capable of 55°C
- 96-well Microplate (500 µL) or desired elution plate
- 2 mL 96-well deep-well plate (Recommend Nunc, Part#278752) or desired plate compatible with the magnetic separation device
- Multichannel pipettes and reagent reservoirs
- Sealing film (Cat# AC1200)
- 100% ethanol
- 100% isopropanol
- Nuclease-free water
- PBS
- Optional: RNase A (10 mg/mL)

Before Starting:

- Prepare SPM Buffer, VHB Buffer, and HDQ Binding Buffer according to the "Preparing Reagents" section on Page 5.
- Set heat block, incubator, or water bath to 70°C.

1. Punch out 3, 5mm blood spots from the filter paper and place at the bottom of 2 mL nuclease-free microcentrifuge tube (not provided).
2. Add 250 µL PBS (not provided) and 20 µL Proteinase K Solution to each sample.
3. Incubate at 55°C in a shaking water bath for 1 hour or until lysis is complete.

Note:

- If a shaking water bath is not available, vortex the samples vigorously every 20-30 minutes. Incomplete lysis may significantly reduce DNA yields.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

4. Add 290 μL AL Buffer. Vortex or pipet up and down 20 times to mix. Proper mixing is crucial for good yield.

Note: For automated protocols tip mixing yields best results and is recommended.

Optional: Add 5 μL RNase A to each sample. Vortex to mix. Let sit at room temperature for 2 minutes.

5. Centrifuge at $\geq 13,000g$ for 5 minutes.
6. Carefully transfer the supernatant to a new 96-well deep-well plate (not provided) without disturbing the filter paper.
7. Add 400 μL HDQ Binding Buffer and 20 μL Mag-Bind® Particles HDQ. Vortex for 10 minutes to mix.

Note:

- HDQ Binding Buffer must be diluted with 100% isopropanol prior to use. Please see Page 5 for instructions. HDQ Binding Buffer and Mag-Bind® Particles HDQ can be prepared as a mastermix. Prepare only what is needed for each run.
- If constant vortexing for 10 minutes is not possible, vortex for 30 seconds every 2 minutes for 10 minutes

8. Place the plate on a magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
9. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
10. Remove the plate from the magnetic separation device.
11. Add 600 μL VHB Buffer.

Note: VHB Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

12. Vortex for 2 minutes.

Note: Complete resuspension of the Mag-Bind® Particles HDQ is critical for obtaining good purity.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

13. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
 14. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
 15. Remove the plate from the magnetic separation device.
 16. Repeat Steps 11-15 for a second VHB Buffer step.
 17. Add 600 µL SPM Buffer.
Note: SPM Buffer must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
 18. Vortex for 2 minutes.
 19. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
 20. Aspirate and discard the cleared supernatant. Do not disturb the Mag-Bind® Particles HDQ.
 21. Select one of the following ethanol removal steps:
 - A. Leave the plate on the magnetic separation device. Add 500 µL nuclease-free water (not provided), leave on magnet for 20-30 seconds, and then aspirate. Do not leave nuclease-free water on Mag-Bind® Particles HDQ for more than 60 seconds. Continue to Step 22.
- OR**
- B. Leave the plate on the magnetic separation device. Wait 1 minute. Remove residual liquid with a pipettor. Dry the Mag-Bind® Particles HDQ for an additional 10 minutes. Continue to Step 22.

Mag-Bind® Blood & Tissue DNA HDQ 96 Kit

22. Remove the plate from the magnetic separation device.
23. Add 50-200 μL Elution Buffer or nuclease-free water to elute DNA from the Mag-Bind® Particles HDQ.

Note: Heat Elution Buffer or nuclease-free water to 70°C to improve yield.
24. Vortex for 5 minutes.

Note: If constant vortexing for 5 minutes is not possible, vortex for 15 seconds every 1-2 minutes for 5 minutes.
25. Place the plate on the magnetic separation device to magnetize the Mag-Bind® Particles HDQ. Let sit at room temperature until the Mag-Bind® Particles HDQ are completely cleared from solution.
26. Transfer the cleared supernatant containing purified DNA to a 96-well microplate (not provided).
27. Store DNA at -20°C.

Note: Blood spots from finger pricks usually contain no more than 50 μL blood, and yield approximately 500 ng to 1 μg DNA. This is usually sufficient for PCR analysis. To obtain higher DNA concentrations, elute with 50 μL pre-heated elution buffer (volumes lower than 50 μL greatly reduce yields).

Troubleshooting Guide

Please use this guide to troubleshoot any problems that may arise. For further assistance, please contact the technical support staff, toll free, at (1-800-832-8896).


Problem	Cause	Solution
Low DNA yield	Incomplete resuspension of Mag-Bind® Particles HDQ	Resuspend the Mag-Bind® Particles HDQ by vortexing vigorously before use.
	Frozen blood samples not mixed properly after thawing	Thaw the frozen blood at room temperature and mix the blood by gently inverting the tube.
	Inefficient cell lysis due to inefficient mixing of AL Buffer/Proteinase K Solution mastermix and sample	Make sure the sample is thoroughly mixed with AL Buffer/Proteinase K Solution master mix within 4 hours of master mix preparation.
	Inefficient binding	Keep Mag-Bind® Particles HDQ suspended in solution with constant vortexing/mixing
	Loss of Mag-Bind® Particles HDQ during operation	Avoid disturbing the Mag-Bind® Particles HDQ during aspiration.
	DNA remains bound to Mag-Bind® Particles HDQ	Increase elution volume and incubation time to 15 minutes; pipet up and down 50 to 100 times.
	DNA washed off	Dilute SPM Buffer by adding appropriate volume of ethanol prior to use (see Page 5 for instructions).
	Ethanol is not added to VHB buffer	Make sure to add ethanol to the VHB Buffer (see Page 5 for instructions).

Troubleshooting Guide

Problem	Cause	Solution
Mag-Bind® Particles HDQ do not completely clear from solution	Too short of magnetizing time	Increase collection time on the magnet.
Gel-like material in the eluted DNA	Blood is too old	Remove the gel-like material by centrifugation; recommend using fresh blood.
		Use 8 mM NaOH as elution buffer.
Problems in downstream applications	Salt carryover	SPM Buffer must be at room temperature.
	Ethanol carryover	Dry the Mag-Bind® Particles HDQ before elution.
	Insufficient washing	Keep Mag-Bind® Particles HDQ suspended during washes with constant vortexing/mixing















Contact Information

To reorder supplies, report a device failure, or complaint, please contact:

 The logo for Omega Bio-tek, Inc. is a stylized black silhouette of a factory or industrial building with three peaks of varying heights.	<p>Manufacturer Omega Bio-tek, Inc. 400 Pinnacle Way Suite 450 Norcross, GA 30071 Website: www.omegabiotek.com Email: info@omegabiotek.com SRN: US-MF-000024148</p>
--	---

Symbols

The following symbols may appear in the instructions for use or on the packaging and labeling:

Picture	Description
 YYYY-MM	Use-by date
	Check components for storage conditions
	Lot number
	Manufacturer
	No additional hazards or not classified as hazardous according to GHS. Also see hazardous symbols as defined in the Precautions Section.
	Website
	Telephone
	Fax
	Email
	LinkedIn
	X
	Facebook
	Recycling Guidance at https://omegabiotek.com/company/recycling
	

Document Revision History

Revision	Description
December 2025 v9.0	<p>Protocol modifications have been made to clarify vortex and incubation steps.</p> <p>Manual has been updated to a new template.</p> <p>Addition of Dried Blood Spot Protocol.</p> <p>Updated recommendations for the magnetic separation device based on the process step.</p>
January 2024 v8.2	<p>Warnings and Safety and Handling Information sections have been added to the manual.</p> <p>A new 24 x 96 prep kit has been added and is now available for purchase.</p> <p>An important statement has been included clarifying how the actual number of preparations is dependent on various factors and may be lower than the number of preparations specified with the kit.</p>
May 2021 v8.1	<p>REACH statement has been added for European Union customers.</p>
March 2020 v8.0	<p>Protocol modifications have been made to the Saliva Protocol to accommodate 500 µL sample volume.</p>
June 2019, v7.0	<p>SPM Wash Buffer has been renamed SPM Buffer. This is a name change only. The formulation has not changed.</p>
January 2019, v6.1	<p>Manual has been updated with a new logo.</p>
August 2018, v6.0	<p>An optional method to dry Mag-Bind® Particles HDQ has been added to Blood, Tissue, Cultured Cells, and Tail Snips extraction protocols.</p> <p>Addition of recommended master mix of AL Buffer and Proteinase K Solution has been added to Blood, Cultured Cells, Saliva, and Buccal Swab extraction protocols.</p>
June 2016, v5.0	<p>M6399 Mag-Bind® Blood DNA HDQ 96 Kit is now Mag-Bind® Blood & Tissue DNA HDQ 96 Kit.</p> <p>Five additional protocols for Tissue, Cultured Cells, Saliva, Buccal Swabs, and Mouse Tail Snips have been included.</p> <p>TL Buffer has been added to the kit for the Tissue Protocol.</p> <p>The volume of Elution Buffer has been increase.</p>

Notices and Disclaimers

For European Union Use.

AL Buffer contains Triton X-100, 2-[4-(2,4,4-trimethylpentan-2-yl)phenoxy]ethanol (CAS 9002-93-1), a substance included in the European Authorisation list (Annex XIV) of REACH Regulation (EC) No 1907/2006. Substances and mixtures used for the purpose of Scientific Research and Development (SR&D) are exempt from authorization requirements if used below 1 tonne per year in volume.

Scientific Research and Development includes experimental research or analytical activities at a laboratory scale such as synthesis and testing of applications of chemicals, release tests, etc. as well as the use of the substance in monitoring and routine quality control or *in vitro* diagnostics.

Trademarks and Licenses

HiBind®, E.Z.N.A.®, MicroElute®, Mag-Bind®, Mag-Binder®, and MB Fit24™ are registered trademarks of Omega Bio-tek, Inc. Qiagen®, QIAvac®, and Vacman® are all trademarks of their respective companies. PCR is a patented process of Hoffman-La Roche. Use of the PCR process requires a license.

Notes:

Notes:

Notes:

For more purification solutions, visit www.omegabiotek.com

AVAILABLE FORMATS



Spin Columns



96-Well Silica Plates



Mag Beads

SAMPLE TYPES



Blood / Plasma



Plasmid



Cultured Cells



Plant & Soil



NGS Clean Up



Tissue




FFPE




Fecal Matter




innovations in nucleic acid isolation

 Omega Bio-tek, Inc.
400 Pinnacle Way, Suite 450
Norcross, GA 30071

 www.omegabiotek.com

 770-931-8400

 770-931-0230

 info@omegabiotek.com

 [omega-bio-tek](https://www.linkedin.com/company/omega-bio-tek)

 [omegabiotek](https://twitter.com/omegabiotek)

 [omegabiotek](https://www.facebook.com/omegabiotek)